Operation Manual

Our eye protection corresponds to the safety requirements of European Regulation (EU) 2016/425 regarding personal safety equipment and was produced according to the standards EN166 / EN170 / EN 171 / EN 172 / EN1731.

Applications:

According to the reference standard. The CE mark and the number of the EN standard are required with the various symbols of the designated area.

Increased strength and symbols for high speed particles	S	Increased strength (12 m/s)
(glass and frame marking)	F	Push with lower energy (45 m/s)
	B	Push with moderate energy (120 m/s)

If the S, F and B symbols on glasses and frames are not the same, the lower of the symbols as the factor for the entire eye protection must be used. The letter T immediately after the strength symbol allows the use of high-energy particles with extremely high temperatures. If the T designation is missing, ambient temperature is assumed.

Symbols for the optical class: (glasses definition)	1	Continuous use
Application icons:	2	Moderate use
	3	Drops or splashes

Standards for application types: (glasses definition)

	Code	Classification
EN 166 (eye protection)	-	-
EN 170 (ultraviolet filter)	2*	1.2 to 5
EN 171 (infrared filter)	4*	1.2 to 10
EN 172 (sun protection filter without infrared specification)	5	1.1 to 4.1
EN 172 (sun protection filter with infrared specification)	6	1.1 to 4.1
EN 1731 (eye/face protection)	-	-

PLEASE READ CAREFULLY THE FOLLOWING HINTS:

CLEANING AND DISINFECTION: The product must be cleaned or disinfected with a cleaning product such as soapy water or a very low-alcohol cleaning solution.

MAINTENANCE: Like all other products for protection, the service life depends on the type of use, treatment and maintenance. If the frame is damaged, replace the entire item. We recommend that glasses and goggles are kept in a suitable case when not in use. In any case, the product must be replaced after a period of 2 years.

WARNING: Glasses with notches and scratched or damaged lenses must be replaced with genuine ones. In certain circumstances, persons with extremely sensitive skin may suffer from allergies caused by material that comes in contact with the skin. A protector worn over prescription eyewear may transmit a shock with a potential risk.

REPLACING GLASSES: Make sure the lenses are the same as the original ones.

TRANSPORT: Eye protection should be transported in the original packaging or an appropriate case.

STORAGE: between 5 and 40°C; humidity < 90%

APPROVALS: Standards and approvals: These products meet the requirements of the applicable standard. Check the product marking for further information on intended use. These products are type approved by independent testing institutes of the European Union and are CE marked to the requirements of Community Directive 89/686/EEC or European Regulation (EU) 2016/425. The applicable legislation can be determined by viewing the Certificate and Declaration of Conformity at www.b-safety24.com.

B-SAFETY GmbH
Oststraße 85, 22844 Norderstedt, Germany
Tel: +49 (0)40-538092-70, Fax: +49 (0)40-538092-86
Web: www.b-safety24.com, E-mail: info@b-safety24.com

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